

# ORBITAL SHAKER ATTACHMENT ACCESSORY FOR THE ENVIRO-GENIE® / INCUBATOR-GENIE®

Installation and Operating Instructions  
Model SI-1275



1. Place the Orbital Shaker Attachment inside the Enviro-Genie® or Incubator-Genie® chamber, aligning the hex shaft with the hex on the pulley of the Orbital Shaker Attachment. Make sure the two feet are in contact with the chamber base.
2. Use the Enviro-Genie®/Incubator-Genie® Rock Mode for controlling the speed of your Orbital Shaker Attachment. Speed range is 10-100 RPM. Refer to the Enviro-Genie® / Incubator-Genie® manual for Rock Mode operation.
3. Place all loads centered on Orbital Shaker platform. Please note, platform capacity is 2 Kg (4.4lb).

 **CAUTION!** Remove from Enviro-Genie® / Incubator-Genie® before cleaning. Do not immerse.

 **CAUTION!** Do not use the Orbital Shaker Attachment in applications such as mixing of flammable materials or where transfer of mechanical energy to glass apparatus could lead to breakage..

 **CAUTION!** Failure to follow operating instructions can compromise the user's safety.

## CARE AND HANDLING

The Orbital Shaker Attachment should be given the care normally required for any appliance. Avoid wetting or unnecessary exposure to fumes. The finish can be washed with a damp cloth and soap or mild detergents. Keep the unit clean by immediately blotting any spills.



80 Orville Drive, Suite 102, Bohemia, NY 11716 U.S.A.  
Phone: (631) 567-4700 • Fax: (631) 567-5896 Toll Free: 888-850-6208  
Customer Service: [custsvc@scientificindustries.com](mailto:custsvc@scientificindustries.com)  
[www.scientificindustries.com](http://www.scientificindustries.com)